## **HAMAMATSU**

# PHOTOMULTIPLIER TUBE R2949

(Low Dark Counts Type of R928)

## Modified R928, 8 mm × 6 mm Photocathode Size Low Dark Counts, Extended Red Multialkali, High Sensitivity For Photon Counting Applications

The R2949 is a 28mm (1-1/8 inch) diameter, 9-stage, side-on type photomultiplier tube having the same extended red multialkali photocathode as the R928. The R2949 features very low dark counts, extremely high quantum efficiency, high gain, good S/N ratio and wide spectral response from UV to near infrared.

The R2949 is well suited for use in various low-level photometries such as general single photon counting applications and fluorescence life time measurement.

The R2949 is directly interchangeable with the R928.

## **FEATURES**

●Low Dark Counts	300 s <sup>-1</sup> (at 25 °C)
●Low Dark Current 2 nA (after 30 mir	n storage in darkness)
●Wide Spectral Response	. 185 nm to 900 nm
●High Cathode Sensitivity	
Luminous	200 μA/lm
Radiant at 400 nm	68 mA/W
●High Anode Sensitivity (at 1000 V)	
Luminous	2000 A/Im
Radiant at 400 nm	6.8 × 10 <sup>5</sup> A/W
●Low Drift and Hysteresis	

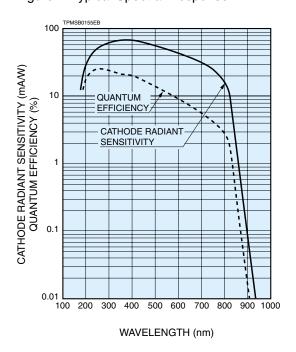


Figure 1: Typical Spectral Response

## SPECIFICATIONS

#### **GENERAL**

Pa	rameter	Description/Value	Unit
Spectral Respo	onse	185 to 900	nm
Wavelength of	Maximum Response	400	nm
Photocathode	Material	Multialkali	_
Priotocatriode	Minimum Effective Area	8×6	mm
Window Materi	al	UV glass	_
	Secondary Emitting Surface	Multialkali	_
Dynode	Structure	Circular-cage	_
	Number of Stages	9	_
Direct Interelectrode	Anode to Last Dynode	4	pF
Capacitances	Anode to All Other Electrodes	6	pF
Base		11-pin base JEDEC No. B11-88	_
Weight		Approx. 45	g
Operating Amb	ient Temperature	-30 to +50	°C
Storage Tempe	rature	-30 to +50	°C
Suitable Socke	t	E678–11A (Sold Separately)	_
Suitable Socke	t Assembly	E717–63 (Sold Separately)	_



Subject to local technical requirements and regulations, availability of products included in this promotional material may vary. Please consult with our sales office.

Information furnished by HAMAMATSU is believed to be reliable. However, no responsibility is assumed for possible inaccuracies or omissions. Specifications are subject to change without notice. No patent rights are granted to any of the circuits described herein. ©2006 Hamamatsu Photonics K.K.

## PHOTOMULTIPLIER TUBES R2949

## **MAXIMUM RATINGS (Absolute Maximum Values)**

	Parameter	Value						
Supply Voltage	Between Anode and Cathode	1250	V					
	Between Anode and Last Dynode	250	V					
Average Anode Curr	rent <sup>A</sup>	0.1	mA					

#### CHARACTERISTICS (at 25 °C)

	Parameter		Min.	Тур.	Max.	Unit
	Quantum Efficiency	at 255 nm	_	25.3	_	%
	Luminous B		140	200	_	μ <b>A</b> /lm
		at 194 nm	_	18	_	mA/W
		at 254 nm	_	52	_	mA/W
Cathode Sensitivity	Radiant	at 400 nm	_	68	_	mA/W
·		at 633 nm	_	41	_	mA/W
		at 852 nm	_	3.5	_	mA/W
	Red/White Ratio C		0.1	0.3	_	_
	Blue Sensitivity Inde	ex <sup>D</sup>	_	7.5	_	_
	Luminous <sup>E</sup>		1000	2000	_	A/lm
		at 194 nm	_	1.8 × 10 <sup>5</sup>	_	A/W
Anada Canaitivity	Radiant	at 254 nm	_	5.2 × 10 <sup>5</sup>	_	A/W
Anode Sensitivity		at 400 nm	_	6.8 × 10 <sup>5</sup>	_	A/W
		at 633 nm	_	4.1 × 10 <sup>5</sup>	_	A/W
		at 852 nm	_	3.5 × 10 <sup>4</sup>	_	A/W
Gain <sup>E</sup>			_	$1.0 \times 10^{7}$	_	_
I Anode Dark Count □		at +25 °C	_	300	500	S <sup>-1</sup>
		at -20 °C	_	3	_	S <sup>-1</sup>
Anode Dark Current F	(After 30 min Storage	in Darkness)	_	2	25	nA
ENI (Equivalent Noise	Input) <sup>G</sup>		_	1.2 × 10 <sup>-16</sup>	_	W
Anoda Pulsa Risa T		īme <sup>H</sup>	_	2.2	_	ns
Time Response <sup>E</sup>	Electron Transit Tim	e <sup>J</sup>	_	22	_	ns
Anada Current Ctability K	Light Hysteresis		_	0.1	_	%
Anode Current Stability K	Voltage Hysteresis		_	1.0	_	%

#### **NOTES**

- A: Averaged over any interval of 30 seconds maximum.
- B: The light source is a tungsten filament lamp operated at a distribution temperature of 2856K. Supply voltage is 100 V between the cathode and all other electrodes connected together as anode.
- C: Red/White ratio is the quotient of the cathode current measured using a red filter (Toshiba R-68) interposed between the light source and the tube by the cathode current measured with the filter removed under the same conditions as Note B.
- D:The value is cathode output current when a blue filter (Corning CS 5-58 polished to 1/2 stock thickness) is interposed between the light source and the tube under the same condition as Note B.
- E: Measured with the same light source as Note B and with the anode-tocathode supply voltage and voltage distribution ratio shown in Table 1.

#### **Table 1:Voltage Distribution Ratio**

Electrode	K	D	y1	/1 Dy		2 Dy3		Dy4		Dy5		Dy6		Dy7		Dy8		Dy9			Р	
Distribution Ratio		1		1	1		1			1		1	1		1		1		-	1		
Supply Voltage: 1000 V,					ı	K: (	Ca	th	od	le,		Dy	: E	) yr	10	de	,	Р	: A	no	de	

- F: Measured with the same supply voltage and voltage distribution ratio as Note E after 30 minute storage in darkness.
- G:ENI is an indication of the photon-limited signal-to-noise ratio. It refers to the amount of light in watts to produce a signal-to-noise ratio of unity in the output of a photomultiplier tube.

$$ENI = \frac{\sqrt{2q \cdot ldb \cdot G \cdot \Delta f}}{S}$$

where  $q = Electronic charge (1.60 \times 10^{-19} coulomb)$ .


ldb = Anode dark current (after 30 minute storage) in amperes.

G = Gain.

 $\Delta f$  = Bandwidth of the system in hertz. 1 hertz is used.

S = Anode radiant sensitivity in amperes per watt at the wavelength of peak response.

- H:The rise time is the time for the output pulse to rise from 10 % to 90 % of the peak amplitude when the entire photocathode is illuminated by a delta function light pulse.
- J: The electron transit time is the interval between the arrival of delta function light pulse at the entrance window of the tube and the time when the anode output reaches the peak amplitude. In measurement, the whole photocathode is illuminated.
- K: Hysteresis is temporary instability in anode current after light and voltage are applied.

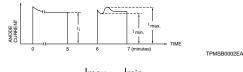

Hysteresis = 
$$\frac{I_{\text{max.}} - I_{\text{min.}}}{I_{\text{i}}} \times 100 \text{ (%)}$$

## HAMAMATSU

#### (1)Light Hysteresis

The tube is operated at 750 V with an anode current of 1  $\mu$ A for 5 minutes. The light is then removed from the tube for a minute. The tube is then re-illuminated by the previous light level for a minute to measure the variation.

Figure 2: Anode Luminous Sensitivity and Gain Characteristics


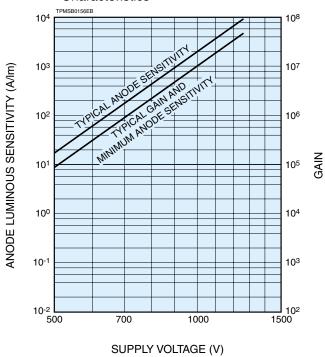


Figure 4: Typical Temperature Coefficient of Anode Sensitivity

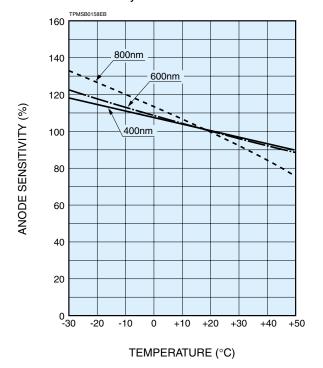

#### (2) Voltage Hysteresis

The tube is operated at 300 V with an anode current of 0.1  $\mu$ A for 5 minutes. The light is then removed from the tube and the supply voltage is quickly increased to 800 V. After a minute, the supply voltage is then reduced to the previous value and the tube is re-illuminated for a minute to measure the variation.

Figure 3: Typical Time Response



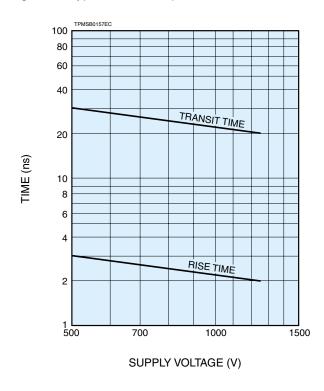


Figure 5: Typical Temperature Characteristic of Dark Counts

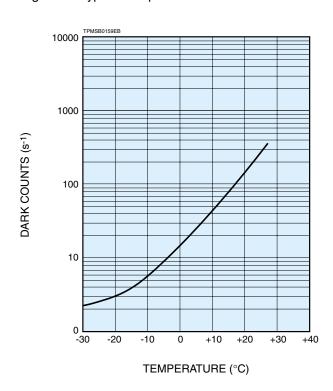


## PHOTOMULTIPLIER TUBES R2949

Figure 6: Single Photon Pulse Height Distribution

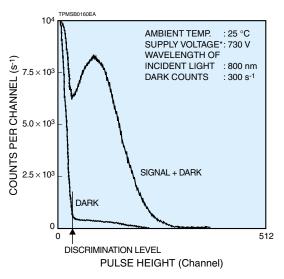

 $<sup>^{\</sup>star}$  The supply voltage is set at gain to be  $1\times10^{6}$ 

Figure 7: Dimensional Outline and Basing Diagram (Unit: mm)

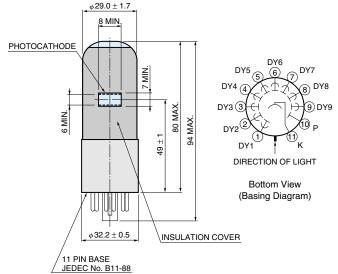


TPMSA0016EA

Figure 7: Accessories (Unit: mm)


Sold Separately


#### Socket E678-11A

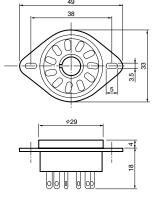

TACCA0064EA

#### D Type Socket Assembly E717-63

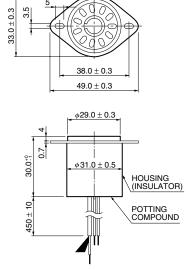

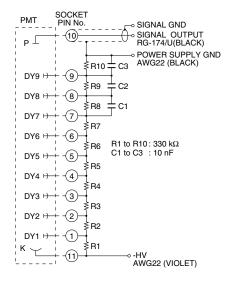

TACCA0002EH

#### Warning-Personal Safety Hazards

Electrical Shock–Operating voltages applied to this device present a shock hazard.

## HAMAMATSU

WEB SITE www.hamamatsu.com

HAMAMATSU PHOTONICS K.K., Electron Tube Division

314-5, Shimokanzo, Iwata City, Shizuoka Pref., 438-0193, Japan, Telephone: (81)539/62-5248, Fax: (81)539/62-2205

U.S.A.: Hamamatsu Corporation: 360 Foothill Road, P. O. Box 6910, Bridgewater. N.J. 08807-0910, U.S.A., Telephone: (1)908-231-0960, Fax: (1)908-231-1218 E-mail: usa@hamamatsu.com
Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany. Telephone: (49)8152-2575-0, Fax: (49)8152-2658 E-mail: info@hamamatsu.de
France: Hamamatsu Photonics France S.A.R.L.: 19, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: (33)1 69 53 71 00, Fax: (33)1 69 53 71 10 E-mail: info@hamamatsu.fr
United Kingdom: Hamamatsu Photonics UK Limited: 2 Howard Court, 10 Tewin Road Welwyn Garden City Hertfordshire AL7 18W, United Kingdom, Telephone: 44-(0)1707-294888, Fax: 44(0)1707-325777 E-mail: info@hamamatsu.co.uk
North Europe: Hamamatsu Photonics Norden AB: Smidesvägen 12, SE-171-41 SOLNA, Sweden, Telephone: (46)8-509-031-00, Fax: (46)8-509-031-01 E-mail: info@hamamatsu.se
Italy: Hamamatsu Photonics Italia: S.R.L.: Strada della Moia, 1/E, 20020 Arese, (Milano), Italy, Telephone: (39)02-935 81 733, Fax: (39)02-935 81 741 E-mail: info@hamamatsu.it

TPMS1045E02
JUL. 2006. IP

<sup>\*</sup> Hamamatsu also provides C4900 series compact high voltage power supplies and C6270 series DP type socket assemblies which incorporate a DC to DC converter type high voltage power supply.